

**FAST, EASY, GENTLE!**

# pluriBead®

**Application Example**

**Isolation of rare cell types:  
Circulating tumor cells**

**pluriSelect  
USA**

Spring Valley, CA 91977  
USA

Phone: 619-928-9265  
support.usa@pluriselect.com  
sales.usa@pluriselect.com

**www.pluriselect.com**


**pluriSelect  
Worldwide**

Deutscher Platz 5c  
04103 Leipzig  
Germany

Phone: +49 341 333858-0  
support@pluriselect.com  
sales@pluriselect.com

V 10

 **pluriSelect**

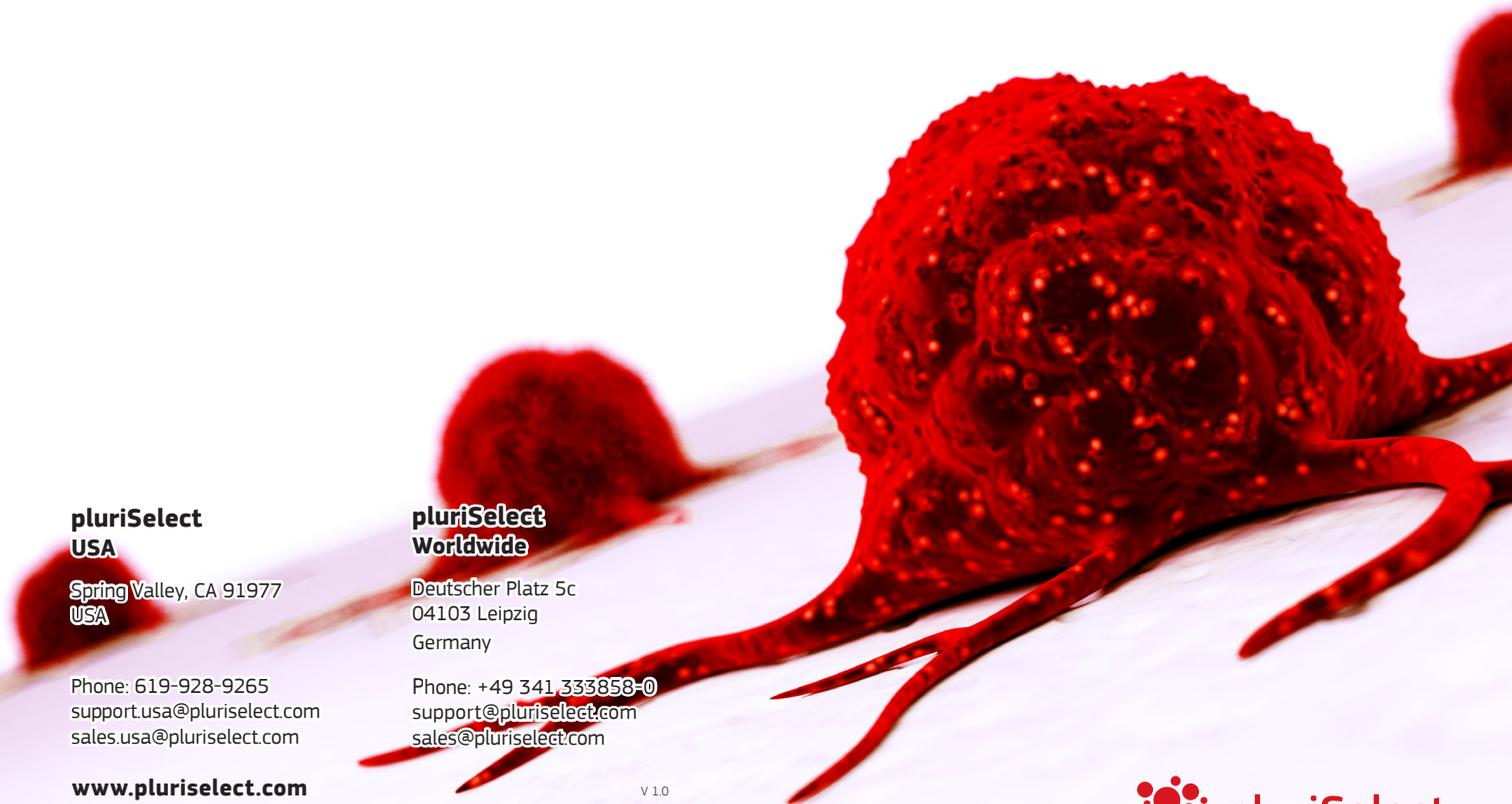


# Isolation of rare cell types: Circulating tumor cells

## Introduction

Here we are presenting a new approach to colon carcinoma circulating tumor cell (CTC) screening using pluriBead® carrying a tumor-associated EpCAM antibody.

This method is based on a non-magnetic cell separation technology. It does not require any sample pre-treatment. EpCAM-pluriBead® can be added directly to a whole blood sample. The method is also suitable for single cell isolation from different biological fluids. Moreover, its sensitivity can be additionally increased by raising the sample volume. The bound EpCAM-positive colon carcinoma cells can be easily involved in further molecular-genetic experiments that aim to detect of their mutation status. In case of colon carcinoma, K-ras mutation status is a predictive tool of response anticancer therapy. Thus, the new method can be considered as a fast and effective instrument for early cancer diagnostics.

## Materials and methods

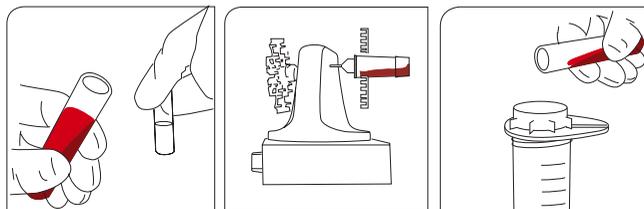

### A: Labeling


Add EpCAM-pluriBead® to your sample.

### B: Incubation

30 minutes of gentle incubation (recommended with pluriPlix®).

### C1: Separation

Captured target cells are separated via appropriate sieves.

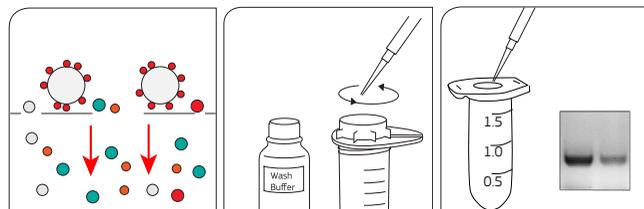

### C2: Separation

Target cells bound on pluriBead® stay on top of the sieve. The rest runs through.

### D: Washing & lysis

Use wash buffer to clean sieve. Lyse cells with Trizol®.

### E: Processing

Approaches for the study of cancer cells:  
- RNA/DNA isolation  
- Cell culture experiments

## Results


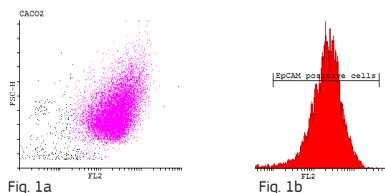


Fig. 1a and 1b: Flow cytometric analysis of Caco-2 cells, staining with anti-EpCAM-PE.

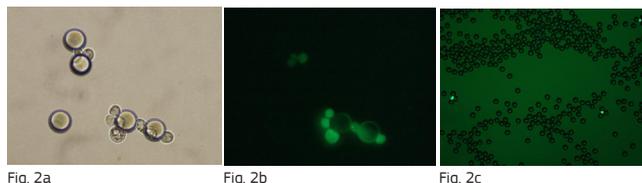

2a) Colon carcinoma Caco-2 cells captured by EpCAM-pluriBead®.

2b) Captured colon carcinoma Caco-2 cells, staining with calcein.

2c) Captured colon carcinoma Caco-2 cells spiked in 30 ml whole blood with 0.5 million EpCAM-pluriBead®.

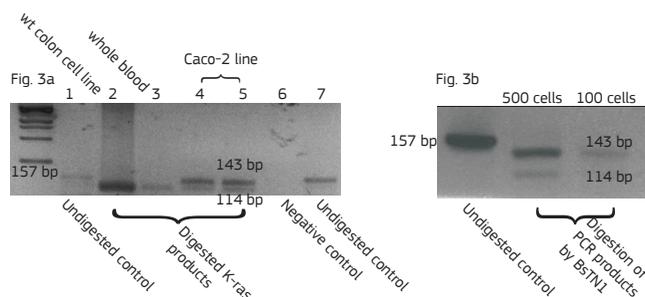


Fig. 3: Screening of colon carcinoma lines for the presence of mutation in 12h codone of K-ras oncogene by PCR using primers that introduce BstNI restriction enzyme sites into PCR products.

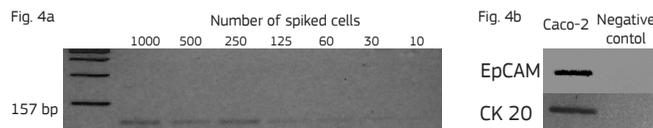


Fig. 4a: Sensitivity of PCR method for the detection of K-ras protooncogene in known number of spiking Caco-2 cells. Fig. 4b: Products of RT-PCR for EpCAM and CK 20 mRNA isolated from the Caco-2 cells, captured by EpCAM-pluriBead® (as a negative control, mRNA from whole blood was used).

## Conclusions and perspectives

Here we developed a new approach for the detection of circulating colon tumor cells in blood using EpCAM-pluriBead®. Captured cells are suitable for further molecular-genetic screening of specific markers, connecting with tumor formation. The developed "in situ immunobeads pcr" method does not require preliminary RNA/DNA isolation and can effectively save the time of analysis. The further aim of this work is to increase the sensitivity of method raising the sample volume, varying the cell tumor lines and/or antibody specificity.