


TwinSpin

Fill – Spin -Drop

Due to movement during transport, the inner tube might be empty. To make sure that there is always Density Gradient Medium inside the tube, perform a quick rotating movement, holding the tube in the upright position. This will fill the inner tube (see figure)

TwinSpin

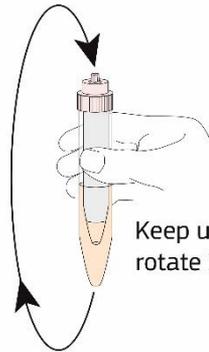

Keep upright,
rotate 1x 360° quickly

TwinSpin was developed for small blood volume (up to 2.50 ml whole blood). TwinSpin for smaller and larger volumes are in the making.

Target cells are concentrated in the opaque ring, more are above and below. Start collecting below.

We also fill TwinSpin with YOUR Density Gradient Medium!!

Pluriselect-USA, Inc. – North America
2030 Gillespie Way 102/103
El Cajon, CA 92020, USA
Phone: +1 619 202-4297
support@pluriselect-usa.com

pluriSelect worldwide
Deutscher Platz 5e
D-04103 Leipzig, Germany
+49 341 333 858 99
info@pluriselect.com